

Special emptying device for grinding bowls *premium line*



Article-description: **Special emptying device for all grinding bowls *premium line* 80 ml, 45 ml, 20 ml volume**
- with 2 sieves – 0.8 + 0.08 mm mesh width -

Article-Number: **50.9900.00**

Description: The emptying device for *premium line* grinding bowls enables a quick and easy separation of grinding balls and suspension after the grinding process.

How it works – look here!



Bowl with balls and suspension after completion of the grinding!



Firmly attach the emptying device with the sieve and seal ring onto the grinding bowl!



Fasten the attachment for the syringe onto the emptying device!



Tense the syringe firmly on the emptying device!



Shake bowl with suspension, draw on the syringe!
(if necessary – pump several times)



The result:
The suspension is in the syringe and the balls are in the bowl!